Human CTP (Cochlin-tomoprotein) Assay Kit - IBL

96 Well

Please read carefully this instruction prior you use this assay kit.

INSTRUCTIONS FOR USE

This product is for research use only and is not intended for diagnostic use.

KIT COMPONENT

1	Precoated plate	96Well x 1
2	Labeled antibody conc. (60X)	0.2mL x 1
3	Standard	1.0mL x 1
4	EIA buffer	30mL x 1
5	Solution for labeled antibody	12mL x 1
6	Chromogen: TMB solution	15mL x 1
7	Stop solution	12mL x 1
8	Wash buffer conc.	50mL x 1

MESUREING SAMPLES

Culture supernatant, Middle Ear Lavage (MEL)

OPERATING PRECATION

- 1 Test samples should be measured soon after collection. For storage of samples, store them frozen and do not repeat freeze/thaw cycles. Thaw the test samples at a low temperature and mix them completely before measurement.
- 2 Test samples should be diluted with "4, EIA buffer" contained in this kit.
- 3 Duplicate measurement of test samples and standards is recommended.
- 4 Standard curve should run for each assay.
- 5 Use test samples in neutral pH range. The contaminations of organic solvent may affect the measurement.
- 6 All reagents should be brought to room temperature (R.T.) and mixed completely and gently before use. After mixing them, make sure of no change in quality of the reagents.
- 7 Use only "8, Wash buffer conc." contained in this kit for washing the precoated plate. Insufficient washing may lead to the failure in measurement.
- 8 Fill the wash buffer each well, invert the plate and make sure the liquid is completely removed by shaking it off if you use a washing bottle. Repeat this washing process several times as instructed in order to avoid any insufficient washing process.
- 9 After remove the wash buffer, tapping the plate against a clean paper towel for completely removing the liquid from the wells and make sure the paper towel is not contact with inside of the wells in this process.
- 10 "6, Chromogen TMB solution" should be stored in the dark due to its sensitivity against light. It should be also avoided contact with metals. Required quantity should be prepared into a collecting container for each use.
- 11 After adding TMB solution into the wells, the liquid in the wells gradually changes the color in blue. In this process the plate should be in dark. Remained TMB solution in the collecting container should not be returned into the original bottle of TMB solution to avoid contamination.
- 12 Measurement of O.D. should be done within 30 minutes after addition of "7, Stop solution".

OPERATION MANUAL AND DOSAGES

- 1. Materials needed but not supplied. Plate reader
 - Test tubes for dilution Deionized water Paper towel Refrigerator

Micropipette and tip Measuring cylinder and beaker Plate washer Collecting container (i.e. clean disposable test tube) Put 230 μ L of standard and gently mix it. Afterword, dilute two fold standard solution in series as a duplicated measurement.

Tube-1 565.0 pg/mL Tube-2 282.5 pg/mL Tube-3 141.3 pg/mL Tube-4 70.6 pg/mL Tube-5 35.3 pg/mL Tube-6 17.7 pg/mL Tube-7 8.8 pg/mL

(4) Preparation of test samples

Dilute test samples with "4, EIA buffer" contained in this kit.*

Conducting pre-assay with different dilutions is recommended for determination of adequate dilution ratio for each testing sample.

3 MEASUREMENT PROCEDURE

(1) Add test sample blank

Determine wells for test sample blank. Put 100µL each of "4, EIA buffer" into the wells.

(2) Add prepared test samples and standard

Put 100 μL prepared test samples and 100 μL prepared standard into appropriate wells.

- (3) Incubation with plate lid (1st reaction).
- (4) Washing
- Wash the plate with the prepared wash buffer and remove all liquid.
- (5) Add prepared labeled antibody
- Put 100 µL prepared labeled antibody into the wells.
- (6) Incubation with plate lid (2nd reaction).
- (7) Washing
- Wash the plate with the prepared wash buffer and remove all liquid completely.
- (8) Add "6, Chromogen TMB solution"

Put 100 µL the TMB solution into the wells.

(9) Incubation in dark

- (10) Add "7, Stop solution"
- Put 100 μL the Stop solution into the wells.
- (11) Determination of optical density (O.D.)

Remove any dirt or drop of water on the bottom of the plate and confirm there is no bubble on the surface of the liquid. Then, measure the both O.D. of standard and the test samples against a test sample blank.

Measurement wavelength: 450 nm. In case of 2 wavelengths:

Main wavelength is 450nm. Sub-wavelength is between 600 and 650 nm.

Table for measurement procedure	Table	for measu	urement r	procedure
---------------------------------	-------	-----------	-----------	-----------

I	able for measurement procedure			
	Test samples	Standard	Test sample blank	
Reagents	Test samples 100 μL	Diluted Standard 100 μL	EIA buffer 100 μL	
1st reaction	Incubation for Overnight 2~8°C with plate lid.			
Washing	4 times (wash buffer more than 350 µL)			
Labeled antibody	100 µL	100 µL		
2nd reaction	Incubation for 30 minutes 2~8°C with plate lid. 5 times (wash buffer more than 350 µL)			
Washing				
TMB solution	100 µL	100 µL	100 µL	
Chromogenic reaction	Incubation for 30 minutes at R.T. (shielded).			
Stop solution	100 µL 100 µL		100 µL	
Measuring O.D.	450 nm / 600~650 nm			

(1) Preparation of wash buffer

Dilute "8, Wash buffer conc." 40 fold with deionized water. The diluted one is used for the assay as a wash buffer. Adjust the required quantities if needed.

(2) Preparation of labeled antibody

Dilute "2, Labeled antibody conc." 60 fold with "5, Solution for labeled antibody" using a prepared collecting container.

(3) Preparation of standard

Prepare 7 test tubes for dilution of the standard and adding 230 μL of the EIA buffer into each tube.

CALCULATION OF TEST RESULT

- 1 Plot the concentration of the standard on the x-axis and its O.D. on the y-axis. Draw a standard curve by applying appropriate regression curve on each plot (4parameter).
- 2 Read the concentration by applying the absorbance of the test samples on a standard curve.
- 3 Calculate the concentration of the test samples by multiplying dilution ratio of test samples on the value.


Manufacturer: Immuno-Biological Laboratories Co., Ltd.

URL: http://www.ibl-japan.co.jp E-mail: do-ibl@ibl-japan.co.jp

Instruction for Use Code No. 27113


Example of standard curve and measured value


Standard (pg/mL)	O.D. (450nm)
565.0	2.577
282.5	1.749
141.3	0.935
70.6	0.481
35.3	0.272
17.7	0.147
8.8	0.082

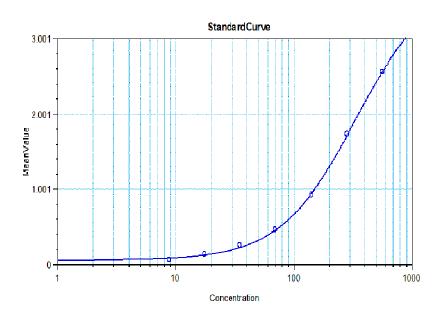

PERFORMANCE CHARACTERISTICS

1. Dilution linearity

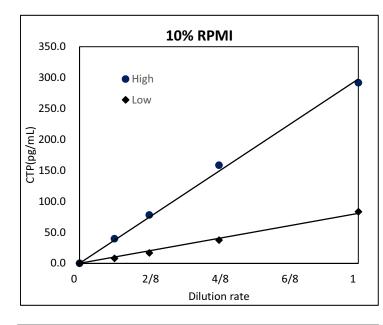

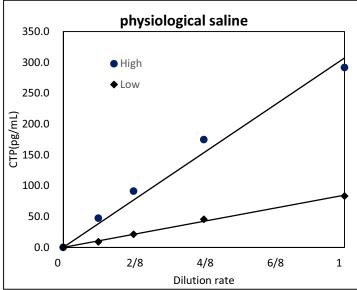

2. Added Recovery Assay

Specimen	Additive Amount (pg/mL)	Theoretical Value (pg/mL)	Measured Value (pg/mL)	%
MEL added recombinant CTP protein	565.00	718.47	711.06	99.0
	282.50	371.05	407.21	109.7
	141.25	184.88	184.55	99.8
	70.63	91.79	97.50	106.2

3. Intra - Assay

Mean Value (pg/mL)	SD (pg/mL)	CV (%)	n
189.32	3.37	1.8	4
43.99	2.54	5.8	4
10.71	0.31	2.9	4

4. Inter - Assay

Mean Value (pg/mL)	SD (pg/mL)	CV (%)	n
179.06	13.12	7.3	12
45.95	3.29	7.2	12
12.51	1.59	12.7	12

PRECAUTION FOR INTENDED USE AND/OR HANDLING

1 Precaution for handling (Hazard prevention)

- (1) Treat the components carefully and wash hands after handling it.
- (2) "7, Stop solution" is a strong acid substance (1N Sulfuric acid). Therefore, it should be careful for the treatment and do not contact your skin and clothes with it. It also needs to pay attention to the disposal of it.

2 Precaution for intended use

- (1) "3, Standard" is lyophilized products. It should be careful to open this vial.
- (2) All reagents should be stored at 2 8°C.
- (3) Precipitation can be seen in "4, EIA buffer", "5, Solution for labeled antibody" and "8, Wash buffer conc.", however, it does not affect its performance.
- (4) Do not mix or replace the reagents with the reagents from a different lot or kit.
- (5) Do not use expired reagents.

3 Precaution for disposal

(1) Dispose used materials after rinsing them with large quantity of water.

STORAGE AND THE TERM OF VALIDITY

Storage Condition: 2 - 8°C The expiry date is specified on the outer box.

PACKAGE UNIT AND PRODUCT NUMBER

Package unit: 96 Well Product number: 27113

CONTACT DETAILS

Immuno-Biological Laboratories Co., Ltd. 1091-1 Naka, Fujioka-Shi, Gunma 375-0005 TEL : 0274-22-2889 FAX : 0274-23-6055

Manufacturer: Immuno-Biological Laboratories Co., Ltd.

URL: http://www.ibl-japan.co.jp E-mail: do-ibl@ibl-japan.co.jp